

MODULE 1: SAMPLE/OBJECTIVE CLEANING

Objective: This module is designed to ensure the proper cleaning of sample and immersion objectives, before and after use. After this module, students should know how to clean in an appropriate and efficient manner.

Step 1)

Remove excess oil from sample/objective by wiping with a **dry** kimwipe.

Step 2)

Place two drops of provided lens cleaner on a **clean** and kimwipe.



Step 3) Begin cleaning the sample/objective using the prepared kimwipe. Start in the center and work your way outwards in a circular motion.

Do this a total of three times:

- 2 times with the kimwipe moistened with lens cleaner and
- 1 time with a dry part of the kimwipe to dry the samples.

Never use the same part of kimwipe during cleaning.

